


Prosonix

Ultrasonic Processing - Prosonitron Technology

Ultrasonic Flow Cells for Laboratory and Plant Scale Processing

30th June 2009

Ultrasonic Processing

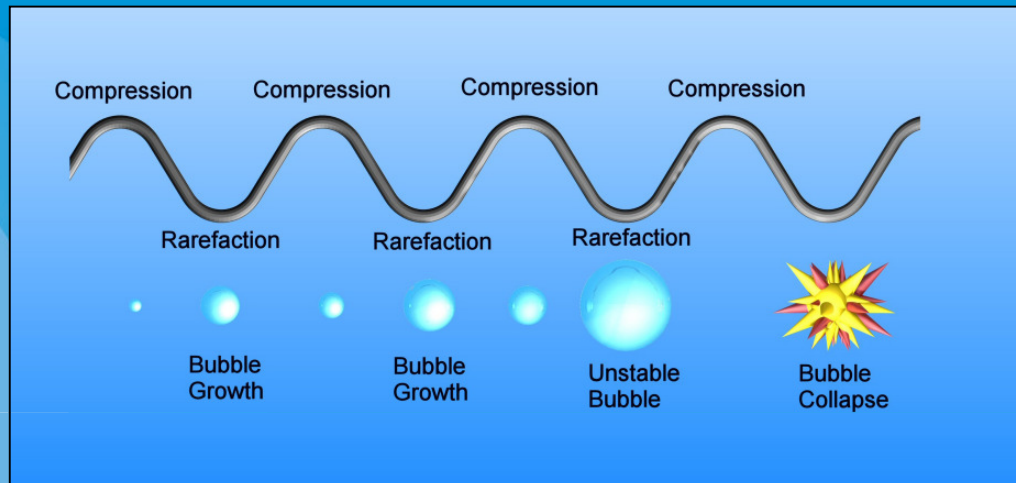


Cavitation Generation with Ultrasound

Sono Nucleation


Sono Milling

Enhanced Mass Transfer

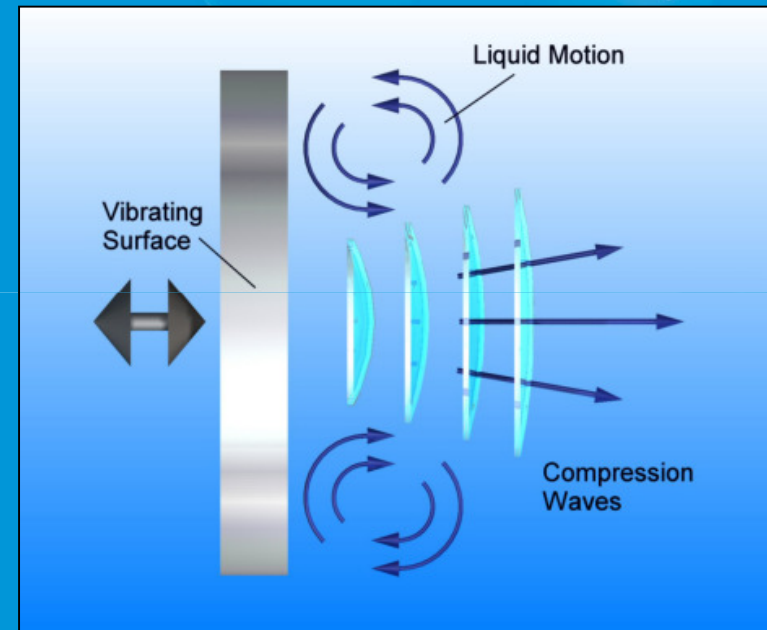


Transmission of Sound Waves

Prosonitron™ Key Feature - Radially Focused Ultrasound

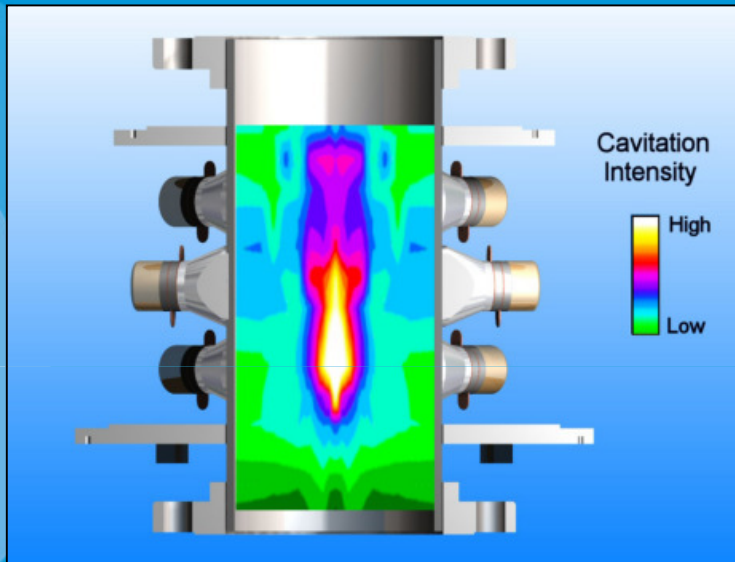

Cavitation Intensity Map of P500 HD
(Taken using probe from NPL)

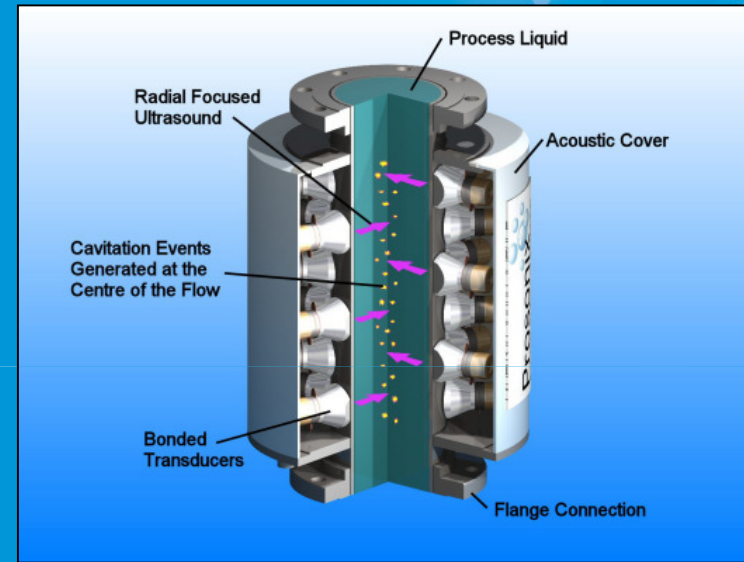


Sound Waves Travel Radially Inward
Building in Intensity to Focus Cavitation at
the Centre of the Cell.



Low surface intensity but high volumetric
intensity allows ultrasonic processing with
minimal surface erosion.

The Prosonitron™ P500 Flow Cell

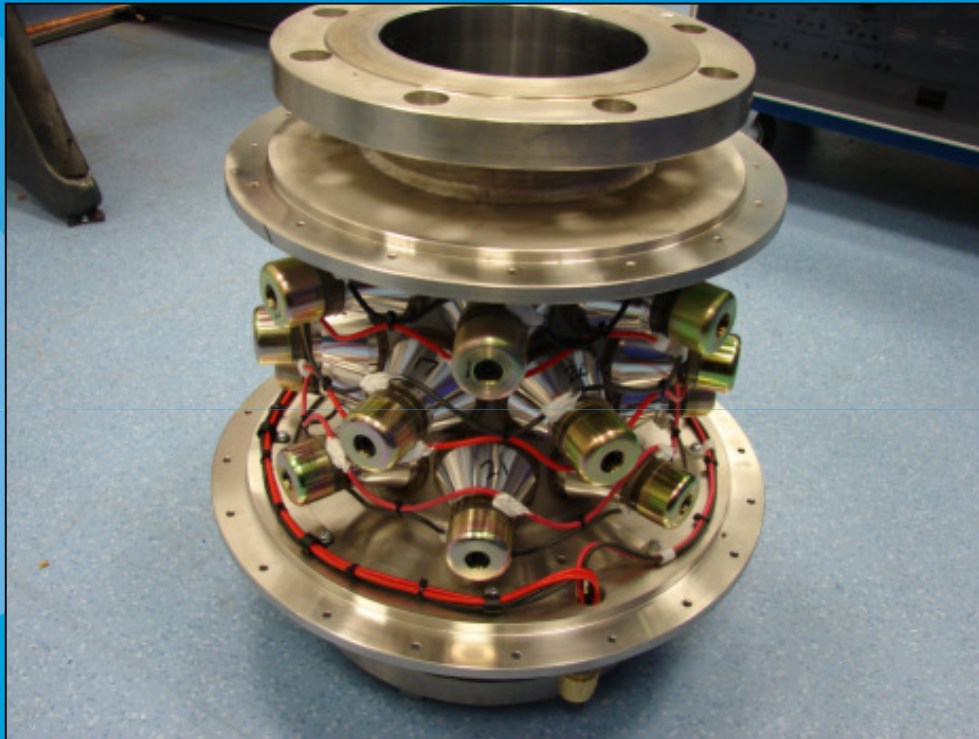

Prosonitron P500 HD flow cell



Generator Cabinet - 2 or 3 Control Cards

Plant Scale Prosonitron Equipment - P500 HD and P750 HD



Prosonitron P750 HD

Prosonitron P500 HD

42 Transducers, 1200 W, 8.2 litre sonicated volume 63 kg

21 Transducers, 600 W, 4.4 litre sonicated volume 49 kg



Plant Installations
Alumina Production

Lab Scale Equipment – SL10 and SL250



Sonolab SL250 – Large Flow Cell

7 Bonded Transducers (50 to 150 W typical) in a similar configuration as the plant scale equipment.

Ideal for 5 to 50 litre batches.



Sonolab SL10 – Mini flow cell

1 Bonded Transducer (5 to 60W typical) delivering power into a 110 ml flow cell volume.

Ideal for 1 to 5 litre batches

Plant Insallations
Alumina Production

Cost Drivers

1. Materials of construction (e.g. Hastelloy C276)
2. Certification of equipment (e.g. ATEX)
3. Aseptic cleaning
4. Size of equipment
 - Configuration
 - Sonication Volume
5. Number of items ordered
6. Commercial terms agreed
 - Warranties
 - Payment terms
 - Service provision

Contact Information

For further information please visit

www.prosonix.co.uk

Or Contact us on

+44 (0) 1865 78 4250